# Delvotest<sup>®</sup> Fast Start III

- EN **Operation Manual**
- DE **Betriebsanleitung**
- FR **Manuel d'utilisation**
- ES **Manual de utilización**
- PT **Operation Manual**
- NL **Gebruikershandleiding**
- TR **Kullanım Kılavuzu**
- IT **Manuale d'uso**
- RU **Руководство по эксплуатации**

EN Although diligent care has been used to ensure that the information provided herein is accurate, nothing contained herein can be construed to imply any representation or warranty for which we assume legal responsibility, including without limitation any warranties as to the accuracy, currency or completeness of this information or of non-infringement of third party intellectual property rights. The content of this document is subject to change without further notice. This document is non-controlled and will not be automatically replaced when changed. Please contact us for the latest version of this document or for further information. Since the user's product formulations, specific use applications and conditions of use are beyond our control, we make no warranty or representation regarding the results which may be obtained by the user. It shall be the responsibility of the user to determine the suitability of our products for the user's specific purposes and the legal status for the user's intended use of our products.

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Version 1.0

### Introduction

Thank you for purchasing our products: Delvotest® Fast Start III. This manual for users contains information and instructions for operation of the incubator. In order to use the instrument properly, please read this manual carefully before use. The Delvotest® Fast Start incubator is designed to be used with Delvotest® Fast range of test for antibiotic residue detection in milk.

### Opening Check

Please check the instrument and adaptor when you first open the packing case. If you find something wrong with the instrument and/or adaptor, contact your local DSM representative.

### Package contents

No.	Item	Type	QTY
1	Incubator	Delvotest® Fast Start	1
2	Adapter	DC12V 5A	1
3	Manual		1

### Safety Precautions

- Read the user manual carefully before operation.
- This product is for indoor use on a levelled surface.
- The metal block heats up during use, do not touch to avoid injuries.
- The operator should not open or repair the instrument. Contact your local DSM representative for assistance.
- The instrument should be used in a location where the temperature is between 5°C (40° F) and 35°C (100° F), it should not be placed near an open window or in direct sunlight. The place should not contain gas or be near a strong disturbing magnetic field. Do not put the instrument in a wet and /or dusty place. The incubator should not be used in a room where antibiotic drugs are stored.
- Power off when you finish your work. Pull out the connector plug when the instrument will not used for long time and prevent dust by storing accordingly.

### Cleaning Instructions

Power off when cleaning incubator. If the instrument gets dirty, clean it with a dry cloth, the metal block should be cleaned regularly with a dry cloth or tissue. Do not use water or any other liquids.

## Operating guide

### Installing the device

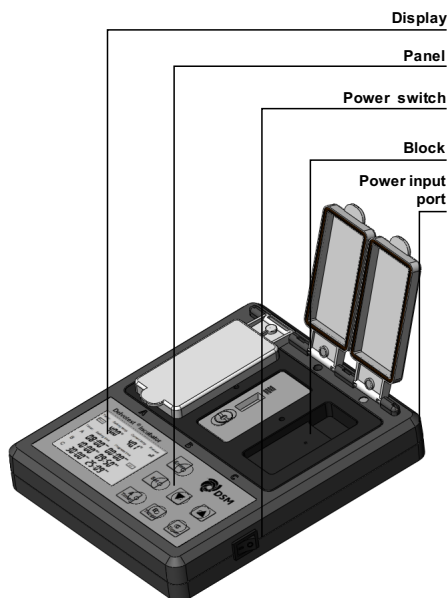
- 1 Place the incubator on a horizontal surface.
- 2 Insert the 12V connector of the adapter into the power connector on the back of the device, and connect the other side of the adapter to the main power supply of 100~240V.
- 3 Turn on the power switch of the device, incubator will start to heat up to the set temperature if preheat function is on. The heating time usually takes 10-15min. See point 4.2.2. pre-heating function.
- 4 Once device reaches to the target temperature, Ready will appear on the top-left of the screen with one beep.

### Keyboard functions

- 1 Set---Press to set the temperature, preheat and time.
- 2 “▲” and “▼”---modify the parameters to set.
- 3 Enter---Save the modified value.
- 4 A Timer / B Timer / C Timer---Short press to start time counting, long press 5 seconds to stop counting. During time counting, press to pause and press again to continue counting.

### Parameter setting

- 1 Press “Set” to enter parameter setting interface. Press “Set” to shift the setting item, press “▲” or “▼” to set parameter value.
- 2 After setting, press “Enter” to save and go back to the main interface. In case of no operation in 10 seconds, device will save the latest value inserted and go back to the main interface automatically.

- In the setting interface, user can set temperature, preheat function and the timer of A, B, C block. Timer of one block can be changed even when other blocks are working.

Note: If the preheat function is "ON", device will automatically heat up to the target temperature after power is switched on. If the preheat function is "OFF", device will only heat up to the target temperature after pressing "Enter".

#### Operation

- After the device is powered on, if "Preheat" is in "ON" status, the incubator will automatically heat to the target temperature. In case "Preheat" is off, press "Enter" to start heating. Once it reaches the target temperature, "Ready" will display on the top left corner of screen with a beep.
- Press the independent buttons "A Timer", "B Timer" and "C Timer" to start time counting for each block.  
Example: Press "A Timer" to start counting for block A. "End" will display with 3 times long beep when time is up. During time counting, press "A Timer" once to pause counting, press again to continue counting. Long press "A Timer" 5 seconds to exit counting.  
Operations on "B Timer" and "C Timer" follows the same method.
- Before the target temperature has been reached, pressing "A Timer" / "B Timer" / "C Timer" will display "Err" with 3 times short beep alarm.
- In case of shut-down or unexpected power-cut, the device will release a short beep alarm.

#### Temperature deviation setting

To be used in case temperature shown on the device is not in line with actual temperature measured with a reference thermometer.

- Press the "Set" and "▼" simultaneously to enter the temperature deviation setting.
- It will display the current deviated value of the temperature on the left of the screen, and display the last-set deviated temperature on the right of the screen.
- Press "▲" and "▼" to set the value of temperature to be deviated and press "Enter" to save and go back to the main interface. The range of temperature deviation allowed is +/- 3°C.
- Long press "Set" 5 seconds to eliminate the deviated value.

Note: If you don't want to set temperature deviation after entering the temperature deviation setting interface, press "Set" and "▼" simultaneously to exit.

#### Troubleshooting

Error	Cause	Solution
No display	No power connection; Power switch failure; Controller failure	Plug in cable on both sides; Check the power switch; Contact vendor
Strong Deviation display temp & actual temp	Sensor broken	Contact vendor
Display "Eoo1" with alarm	Over-heating temperature alarm	Turn off the power and contact vendor
Display "Eoo5" with alarm	Sensor open circuit	Contact vendor
Display "Eoo6" with alarm	Sensor short circuit	Contact vendor
Display "Eoo8" with alarm	Heater failure	Contact vendor
Press invalid	Keyboard failure	Contact vendor

#### Technical data

Model	
Power supply	DC12V 5A
Power capacity	45 W
Average power consumption	10 W
Temperature range	RT+5°C ~ 70°C
Timing range	1sec ~ 99min59 sec
Temperature accuracy	± 2,0°C
Display accuracy	0,1°C
Heating time (from 25°C to 70°C)	≤15min
Ambient temperature	5°C ~ 35°C
Dimensions Device (WxDxH)	188mm x 227mm x 55mm
Dimensions Block (WxDxH)	160mm x 105mm x 18mm
Weight	1,1 kg

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**DSM Food Specialties B.V.**